

DULBECCO'S MINIMAL EAGLE'S MEDIA (DMEM)

13.40 g	DMEM (Gibco 430-2100)
3.70 g	Sodium bicarbonate
2.38 g	Hepes base (Sigma H2275)
0.29 g	Glutamine (Sigma G3126)
0.11 g	Sodium pyruvate (Sigma H2256)
40.0 mg	Gentamycin

Make up to 1 litre with distilled water. Sterilize by filtration (using Millipore's positive pressure system, not by vacuum filtration) and store at 4°C for not more than six weeks.

FETAL CALF SERUM

To heat inactivate calf serum, heat thawed bottles at 56°C for 30 minutes.