

LB (LURIA-BERTANI) MEDIUM

REFERENCE: Molecular Cloning, A Laboratory manual. 1st ed. Maniatis. 1987 p.68.

	<u>HIGH SALT</u>	<u>NORMAL SALT</u>	<u>LOW SALT</u>	<u>NO SALT</u>
Bacto – tryptone	10 g	10 g	10 g	10 g
Bacto – yeast extract	5 g	5 g	5 g	5 g
NaCl	10 g	5 g	0.5 g	-
dH ₂ O	1 litre	1 litre	1 litre	1 litre

For agar plates add 20g agar per litre.
Autoclave all ingredients together to sterilize.

Note: Use low salt or no salt for Pseudomonas